

PUR-SENSE™ Autoclavable and Steam Sterilizable pH Sensors

For additional information, please refer to the Instruction Manuals CD shipped with this product, or visit our website at www.emersonprocess.com/raihome/liquid/.

SPECIFICATIONS

Sensor Type: PUR-SENSE™ 3800, 3800VP, GPHT glass with hemi bulb

Part Number: 3800-01, -02, -03 and 3800VP-01, -02, -03

Range: pH 0 to 14

Temperature Range: up to 105°C (steam up to 140°C) / up to 22 F°
(steam up to 284 F°)

Pressure Range: max 6 bar at 105°C/ max 72 psig (600kPa) at 221 F°)

Temperature sensor: Pt 100

Wetted Materials: Glass, ceramic and EPDM o-ring
USP VI

Reference Junction: Single Ceramic junction

Electrolyte: Gelled polymer

Electrical Connector: Type VP 6 - multipin pH connector
head, suitable for pH and temperature; or coaxial
S8, Pg 13.5 connector

Dimensions: 12 mm OD, shaft length -01= 120 mm ,
-02 = 225 mm , -03 = 325 mm

Remarks: Suitable for all Rosemount pH-instruments. Quality certification includes list of wetted materials and calibration records.



WARNING

Do not exceed temperature and pressure limitations of 140°C (284°F) and 72 psig (600 kPa, 6 bar).

CAUTION

Internal electrolyte fill solution may cause skin or eye irritation.

WARNING



Before removing the sensor, be absolutely certain that the process pressure is reduced to 0 psig and the process temperature is lowered to a safe level!

NOTICE

SENSOR/PROCESS APPLICATION COMPATIBILITY

The wetted sensor materials may not be compatible with process composition and operating conditions. Application compatibility is entirely the responsibility of the user.

STORAGE

1. It is recommended that electrodes be stored in their original shipping containers until needed.
2. Do not store at temperatures below -5°C (23°F).
3. Electrodes should be stored with a protective cap containing KCl solution (PN 9210342).
4. For overnight storage, immerse the sensor in tap water or 4 pH buffer solution.
5. A pH glass electrode does have a limited shelf life of one year.

ELECTRODE PREPARATION

1. Remove electrode from shipping container.
2. Remove the protective boot covering the electrode bulb.
3. Rinse away salt film with clean water, then shake the electrode so that the internal solution fills the bulb, thus removing any air trapped there.



NOTE

Do not allow lubricant to coat electrode bulb or reference junction. If it does, wipe it clean before installation.

INSTALLATION

For sensor orientation, see Figure 2.

For wiring, see see Figures 3 through 8.

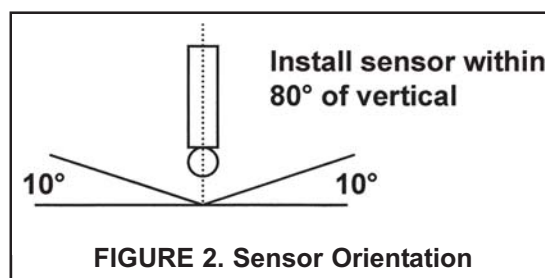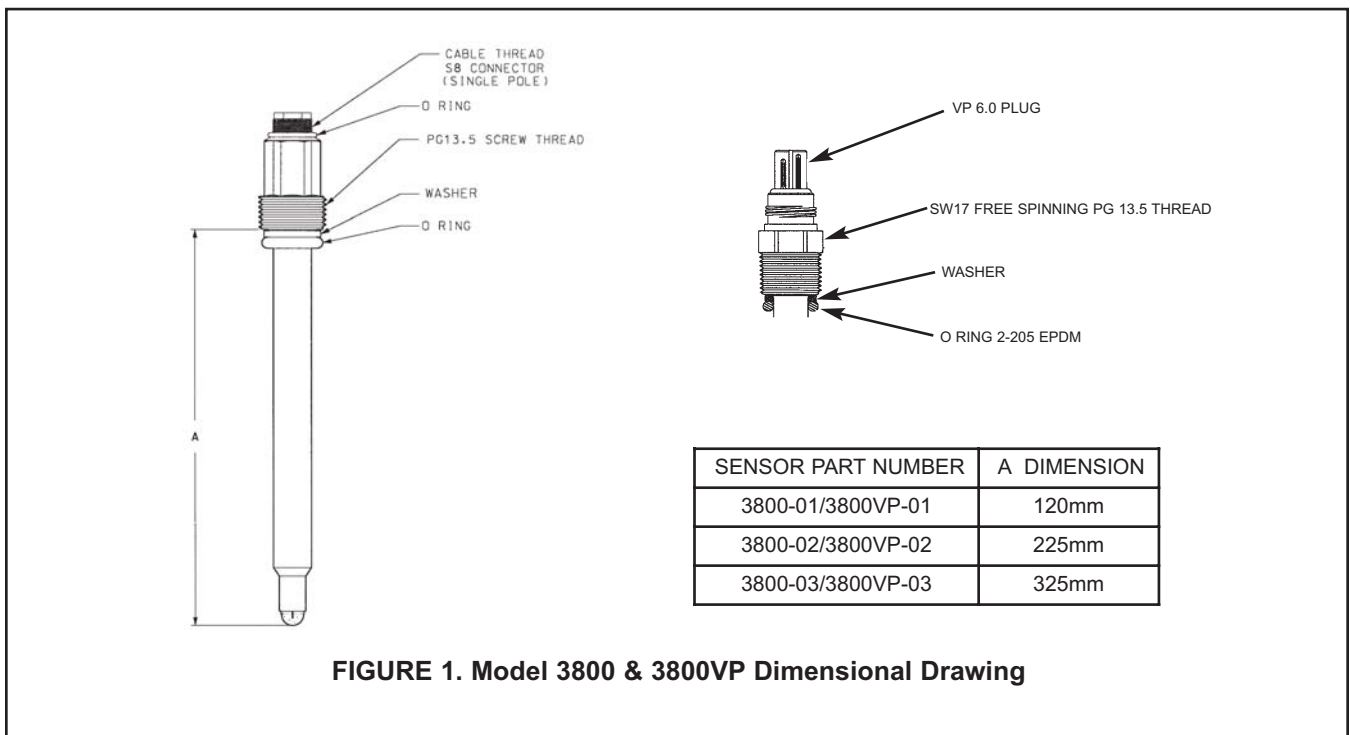

TWO POINT BUFFER CALIBRATION

Select two stable buffer solutions, preferably pH 4.0 and 10.0 (pH buffers other than pH 4.0 and pH 10.0 can be used as long as the pH values are at least two pH units apart).

NOTE

A pH 7 buffer solution reads a mV value of approx. zero, and pH buffers read approximately ± 59.1 mV for each pH unit above or below pH 7. Check the pH buffer manufacturer specifications for millivolt values at various temperatures since it may affect the actual value of the buffer solution mV/pH value.

1. Immerse sensor in the first buffer solution. Allow sensor to equilibrate to the buffer temperature (to avoid errors due to temperature differences between the buffer solution and sensor temperature) and wait for reading to stabilize. Value of buffer can now be acknowledged by analyzer/transmitter.
2. Once the first buffer has been acknowledged by the analyzer/transmitter, rinse the buffer solution off of the sensor with distilled or deionized water.
3. Repeat steps 1 and 2 using the second buffer solution.
4. The theoretical slope value, according to the Nernst equation for calculating pH, is approximately 59.17 mV/pH. Over time the sensor will age, both in the process and in storage, and will result in reduced slope values. To ensure accurate readings, it is recommended that the electrode be replaced when the slope value falls below 47 to 49 mV/pH.

RECOMMENDED pH SENSOR STANDARDIZATION



For maximum accuracy, the sensor can be standardized on-line or with a process grab sample after a buffer calibration has been performed and the sensor has been conditioned to the process. Standardization accounts for the sensor junction potential and other interferences. Standardization will not change the sensor's slope but will simply adjust the analyzer's reading to match that of the known process pH.


MAINTENANCE

Electrodes should respond rapidly. Sluggishness, offsets, and erratic readings are indicators that the electrodes may need cleaning or replacement.



1. To remove oil deposit, clean the electrode with a mild non-abrasive detergent.
2. To remove scale deposits, soak electrodes for 30 to 60 minutes in a 5% hydrochloric acid solution.
3. Temperature effect on life expectancy: If glass electrode life expectancy is 100% @ 25°C (77°F), then it will be approximately 25% @ 80°C (176°F), and approximately 5% @ 120°C (248°F).

WIRING

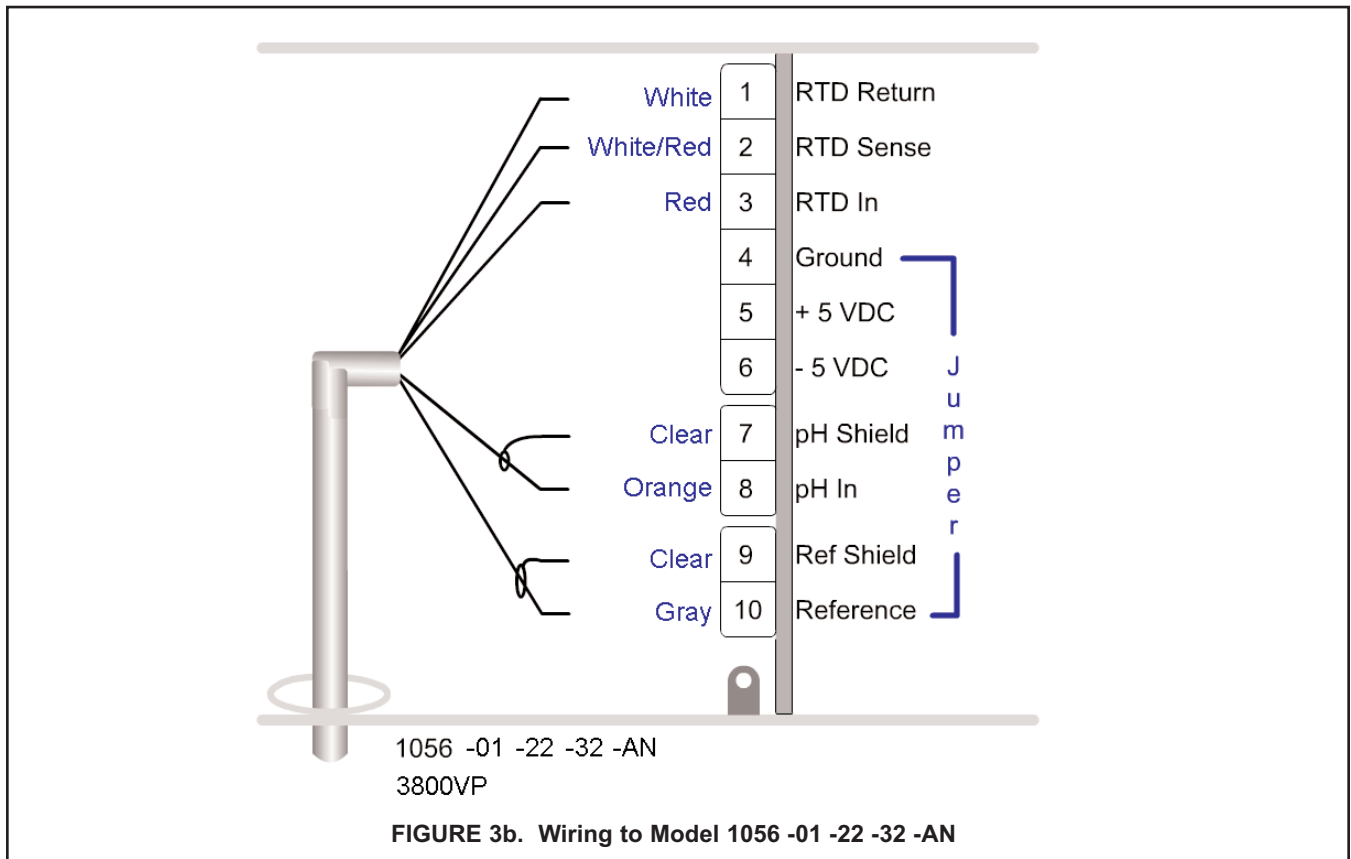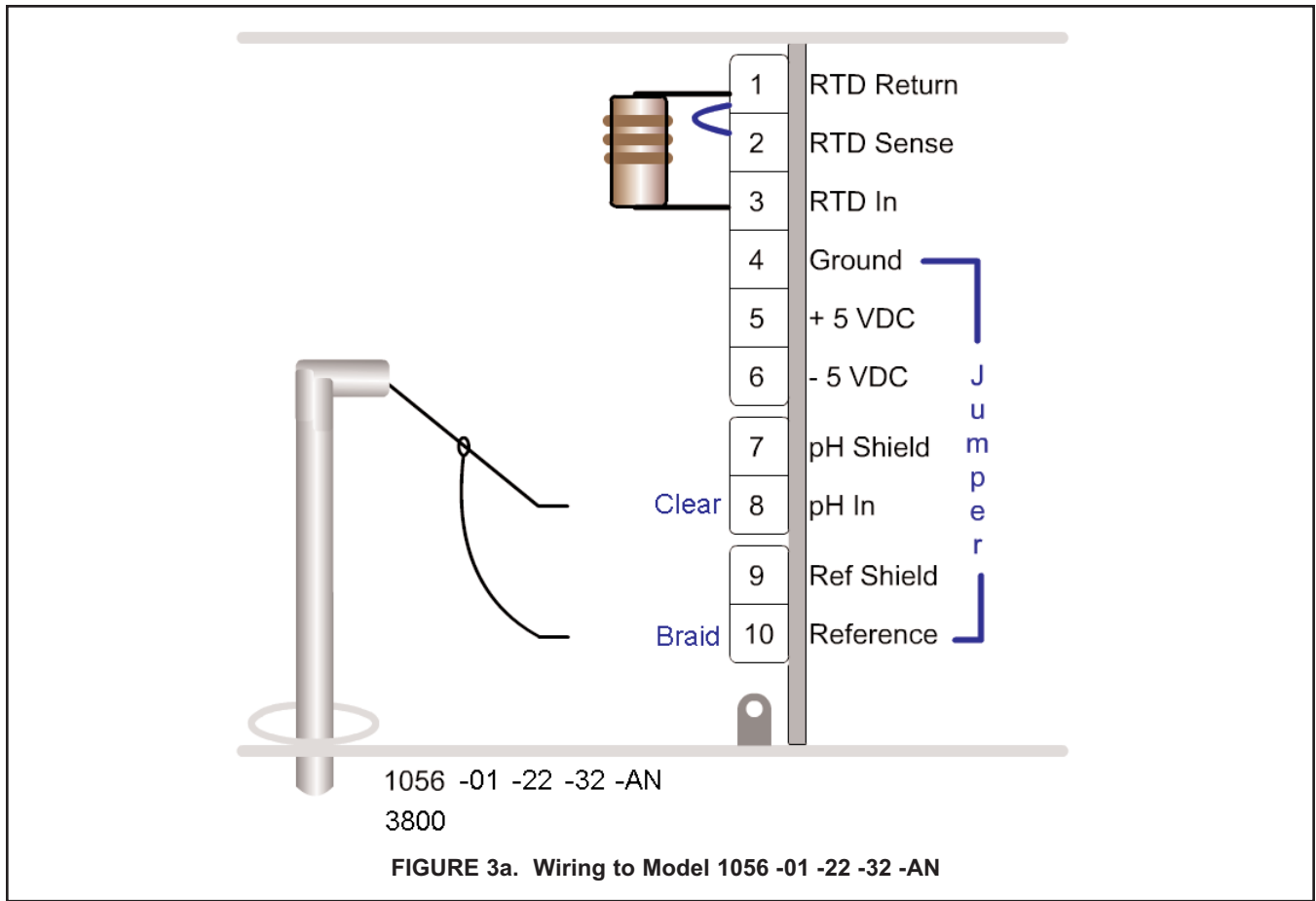

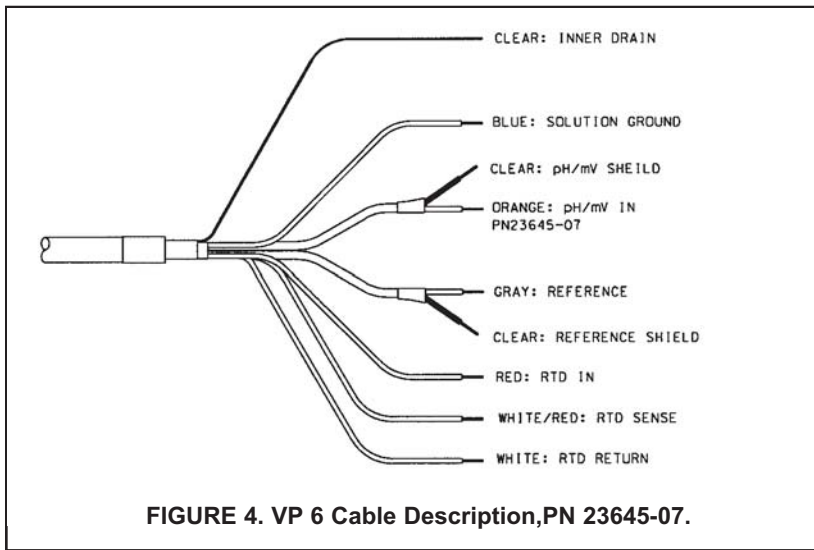

NOTE: For a protective cap is used with holders it must be modified for use with the VP 6.0 cable. The aperture must be enlarged to 7 mm diameter to accommodate this cable.

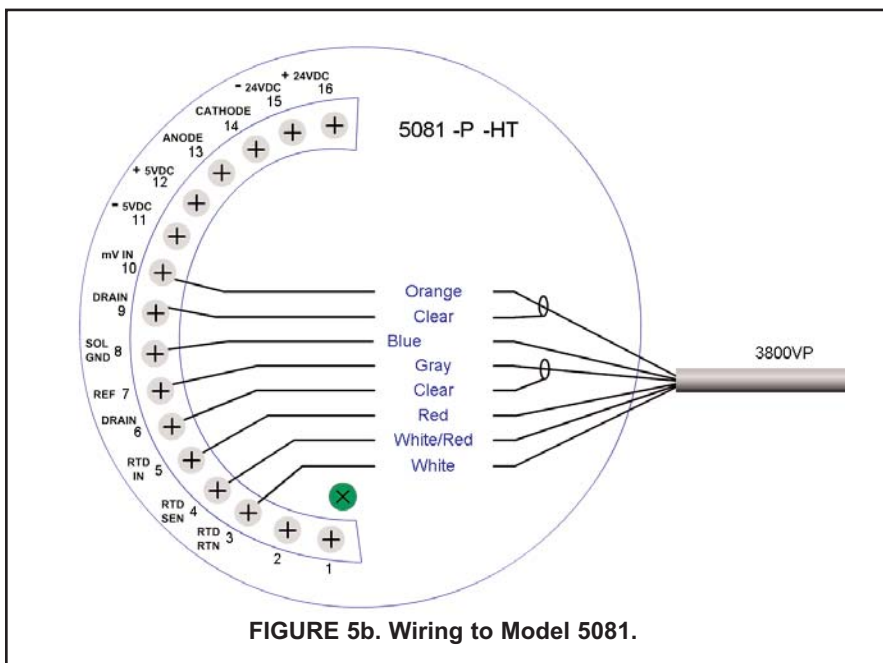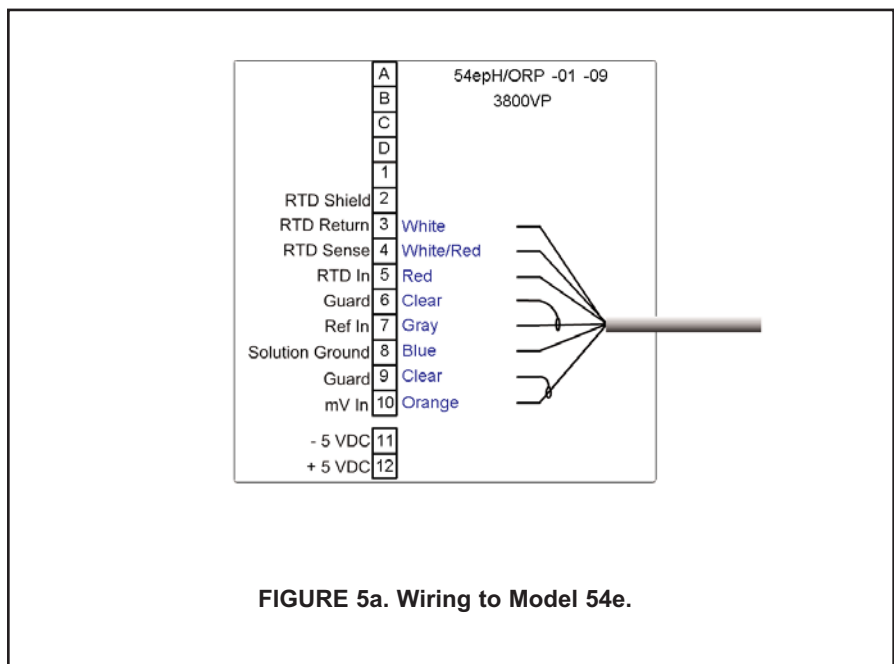

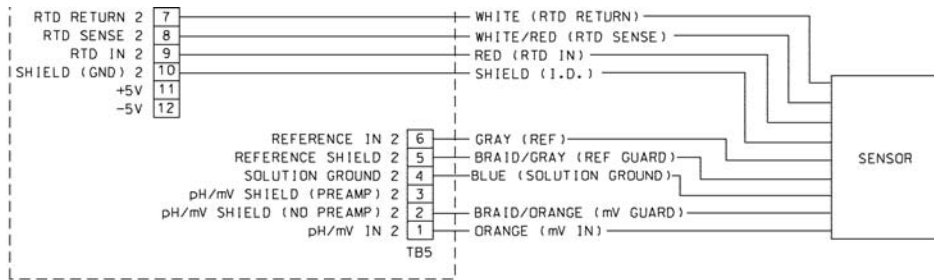


FIGURE 6a. Model 1055-01-10-22 single sensor wiring

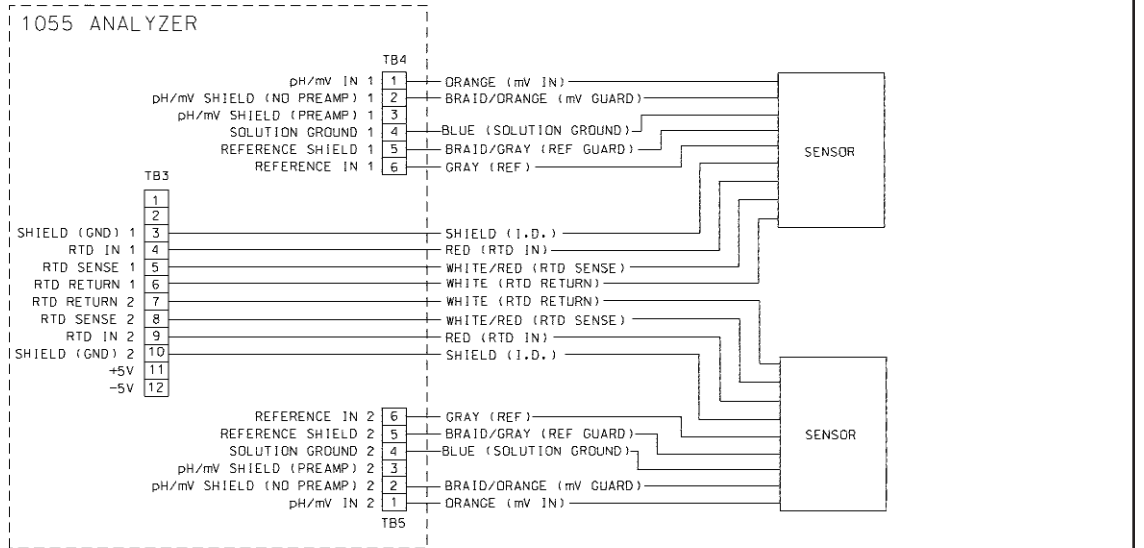


FIGURE 6b. Wiring to Model 1055-01-10-22-32 dual sensor wiring

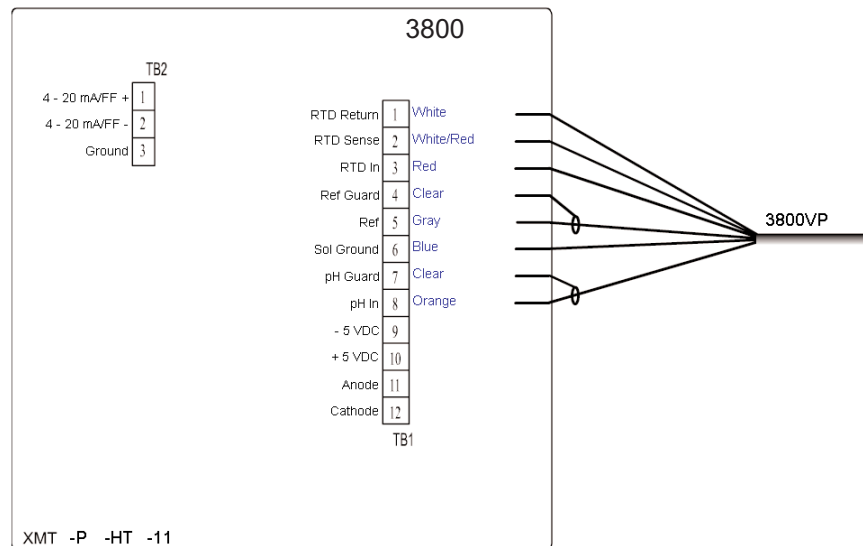


FIGURE 7a. Wiring 3800VP to Model Xmt-P-HT-11 (Pipe Surface Mount)

Note: If no there is no temperature compensation (RTD), then install a 100-ohm resistor between TB1-1 and TB1-3, and install a jumper between TB1-1 and TB1-2.

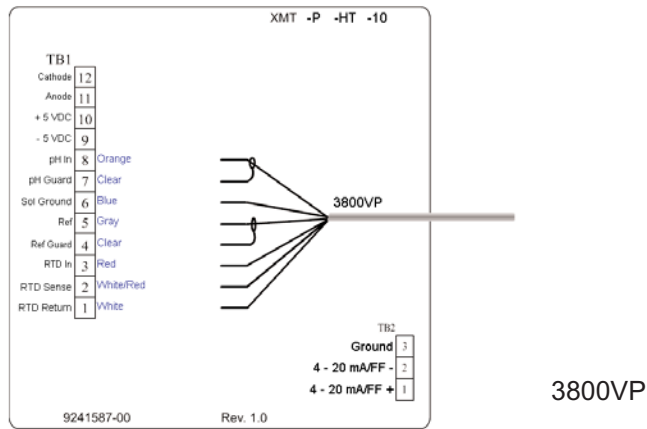


FIGURE 7b. Wiring to Model Xmt-P-HT-10 (Panel Mount)

If no there is no temperature compensation (RTD), then install a 100-ohm resistor between TB1-1 and TB1-3, and install a jumper between TB1-1 and TB1-2.

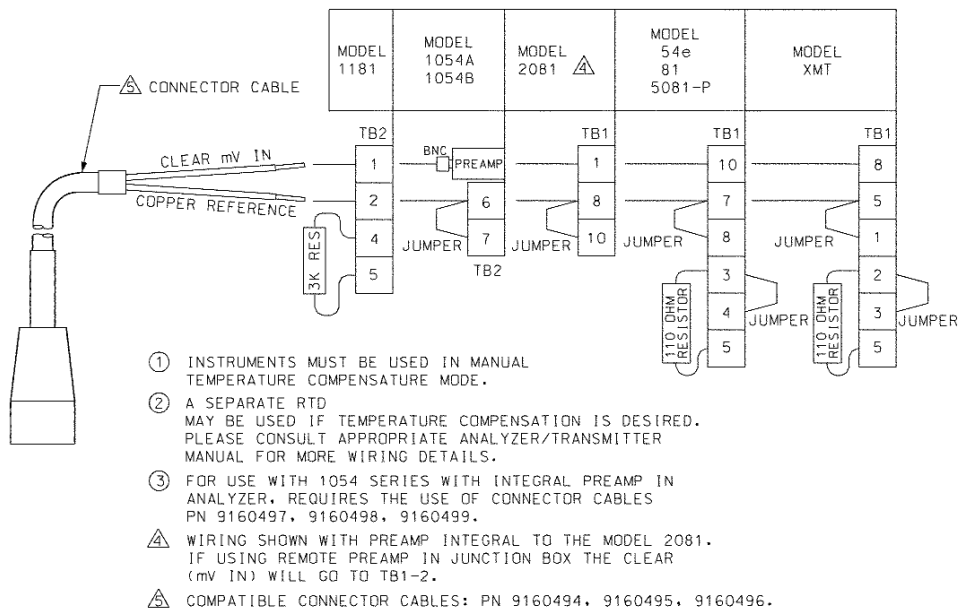


FIGURE 8. Wiring Drawing For Model 3800 with S8 Connector



*The right people,
the right answers,
right now.*

**ROSEMOUNT ANALYTICAL
CUSTOMER SUPPORT CENTER
1-800-854-8257**



Emerson Process Management

Liquid Division

2400 Barranca Parkway
Irvine, CA 92606 USA
Tel: (949) 757-8500
Fax: (949) 474-7250

<http://www.raihome.com>

ON-LINE ORDERING NOW AVAILABLE ON OUR WEB SITE
<http://www.raihome.com>

Specifications subject to change without notice.



Credit Cards for U.S. Purchases Only.

